

Key Features

- Compact unit
- Ideal for microtitre plates and petri dishes
- Laminar flow air circulation
- Digital setting and display of temperature
- Convenient upward opening door
- Stainless steel interior



SI19



SI19

SI19



Incubator, for microtitre plates,

Specially designed for use with microtitre plates, which can be difficult to incubate. In conventional incubators, when 4 or more plates are placed on a shelf, they block the air circulation, which can seriously impair temperature distribution.

The Stuart® SI19 Incubator overcomes this problem with forced air circulation, which blows preheated air across the shelves rather than just through them. Microtitre plates are also prone to drying out during incubation. The SI19 overcomes this with a water tray to increase the humidity within the chamber.

The bright LED display indicates temperature and the encoder control allows rapid and accurate temperature setting. An integral over-temperature protection system automatically tracks the set temperature and controls the heater in the event of a fault. Error codes are displayed if a fault has occurred or if the set temperature has been interrupted.

An acrylic upwards opening door provides excellent access to the interior. The stainless steel chamber features 4 shelves each accommodating 6 plates (or 12 stacked in 2 layers). Shelves are captive and slide out for easy loading.

Technical Specification

| | |
|-------------------------------------|-----------------------|
| Nominal capacity | 20 litres |
| Temperature range | Ambient + 8°C to 80°C |
| Temperature fluctuation (at 37°C) | ±0.5°C |
| Temperature variation | ±0.5°C |
| Display type | LED |
| Display resolution | 0.1°C |
| Internal dimensions, mm (w x d x h) | 250 x 230 x 200 |
| Overall dimensions, mm (w x d x h) | 380 x 380 x 435 |
| Net weight, kg | 24 |
| Electrical supply | 230V, 50Hz, 280W |
| IP Rating | 30 |

Ordering Information

| Model | Description |
|-------|-----------------------------|
| SI19 | Incubator, microtitre plate |

Distribuito da:
Zetalab s.r.l. **Zetalab.it**
Via Umberto Giordano, 5 - 35132 Padova
Tel 049 2021144 - Fax 049 2021143
www.zetalab.it - email: info@zetalab.it